


Electrode to Connector Map

Assembly type	Any 'A' style 16-channel probe on HC16 package.	
Version	V1. Mar 2010	
Connector	Omnetics nano NPD series (A7996-001)	
Mating Connector	Omnetics nano NPD series (A7519-001)	
Compatible Headstages	TDT	RA16CH, LP16CH
	Multi-Channel Systems	Two MPA8I
	Plexon	HST/16o25-18P, HST/16V-G20
	Neuralynx	HS-16
	Alpha Omega	16 mini, TeleSpike
	DataWave Technologies	DW-T-HS16, DW-W-HS16, DW-M-HS16
Compatible Adaptors	TDT	AC-CH
Reference site	Present on limited 16-channel designs. Please check the catalog CAD drawings.	
Note	Electrode sites face up (as shown in the picture).	

CONNECTOR PIN OUT

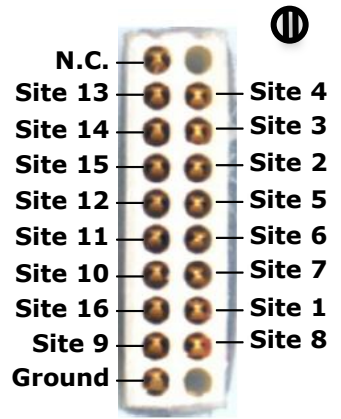

Looking down in the direction as shown by the arrow in the figure below. The inscription may vary on actual probe.

ELECTRODE SITE MAP

