

Radiens Release Notes

Updated June 6, 2024

Radiens Update

A new version of Curate has been released.

<https://www.neuronexus.com/radiens-allego-download-page/>

Curate 1.5.4:

Features/Changes:

- **Protocol Names:** Hitting 'Enter' when typing new protocol name from default name will rename current protocol to the new name, instead of creating a new empty one
When typing in protocols autocomplete field:
 - if "My New Protocol..." is the current selection, it gives an option to save as, which keeps the current protocol graph but saves it under the new name being typed
 - if the current selection is already a custom protocol name, it gives both the Save As option, as well as a Create New option, which creates and opens a new blank protocol
- **Performance:** improvements when loading datasources
- **Bulk Processing:** Bulk Source and Bulk Sink datasource nodes added
- **Status:** Display success or error messages in pop-up upon protocol completion
- **Nodes:** Add context menu on node click with 'Remove node' option
- **Nodes:** 'Open files' button from Data Files tab also creates a Source/Bulk Source node on the protocol, in open Curate tabs
- **Nodes:** Add 'Remove selected nodes' button
- Button styling improvements

Fixes:

- Fix crash when creating new protocol
- Fix crash on reading new generation of sort files
- Improved handling for maximized tabs
- Improvements to conversion from PL2 format
- Improve usability of Remap Sensor node
- Slice Channels node uses site selection state from recording as the initial slicing selection